




β' -Phenyl fentanyl

The Drug Enforcement Administration's Special Testing and Research Laboratory generated this monograph using structurally confirmed reference material.

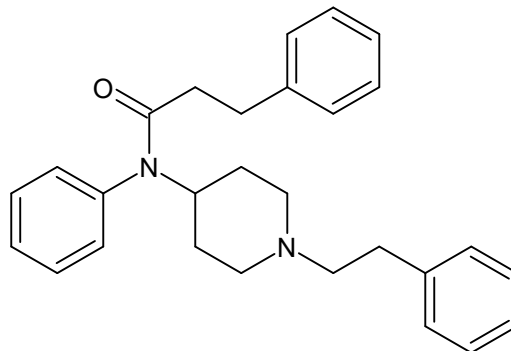

1. GENERAL INFORMATION

IUPAC Name:	<i>N</i> -(1-phenethylpiperidin-4-yl)- <i>N</i> -3-diphenylpropanamide
CAS#:	79279-02-0
Synonyms:	3-phenylpropanoyl fentanyl, hydrocinnamoyl fentanyl, <i>N</i> ,3-diphenyl- <i>N</i> -[1-(2-phenylethyl)piperidin-4-yl]propanamide
Source:	DEA Reference Material Collection
Appearance:	White powder
UV_{max}(nm):	Not determined

2. CHEMICAL AND PHYSICAL DATA

2.1 CHEMICAL DATA

Form	Chemical Formula	Molecular Weight	Melting Point (°C)
Base	C ₂₈ H ₃₂ N ₂ O	412.57	Not Determined
HCl	C ₂₈ H ₃₂ N ₂ O HCl	449.03	Not Determined



β' -Phenyl fentanyl

The Drug Enforcement Administration's Special Testing and Research Laboratory generated this monograph using structurally confirmed reference material.




3. QUALITATIVE DATA

3.1 NUCLEAR MAGNETIC RESONANCE

Sample Preparation: Dilute analyte to ~11 mg/mL in CDCl_3 containing TMS for 0 ppm reference and 1,4-BTMSB- d_4 as quantitative internal standard.

Instrument: 400 MHz NMR spectrometer
Parameters: Spectral width: at least containing -3 ppm through 13 ppm
Pulse angle: 90°
Delay between pulses: 45 seconds

$^1\text{HNMR}$: β' -Phenyl fentanyl; LOT # 0519541-4; CDCl_3 ; 400MHz

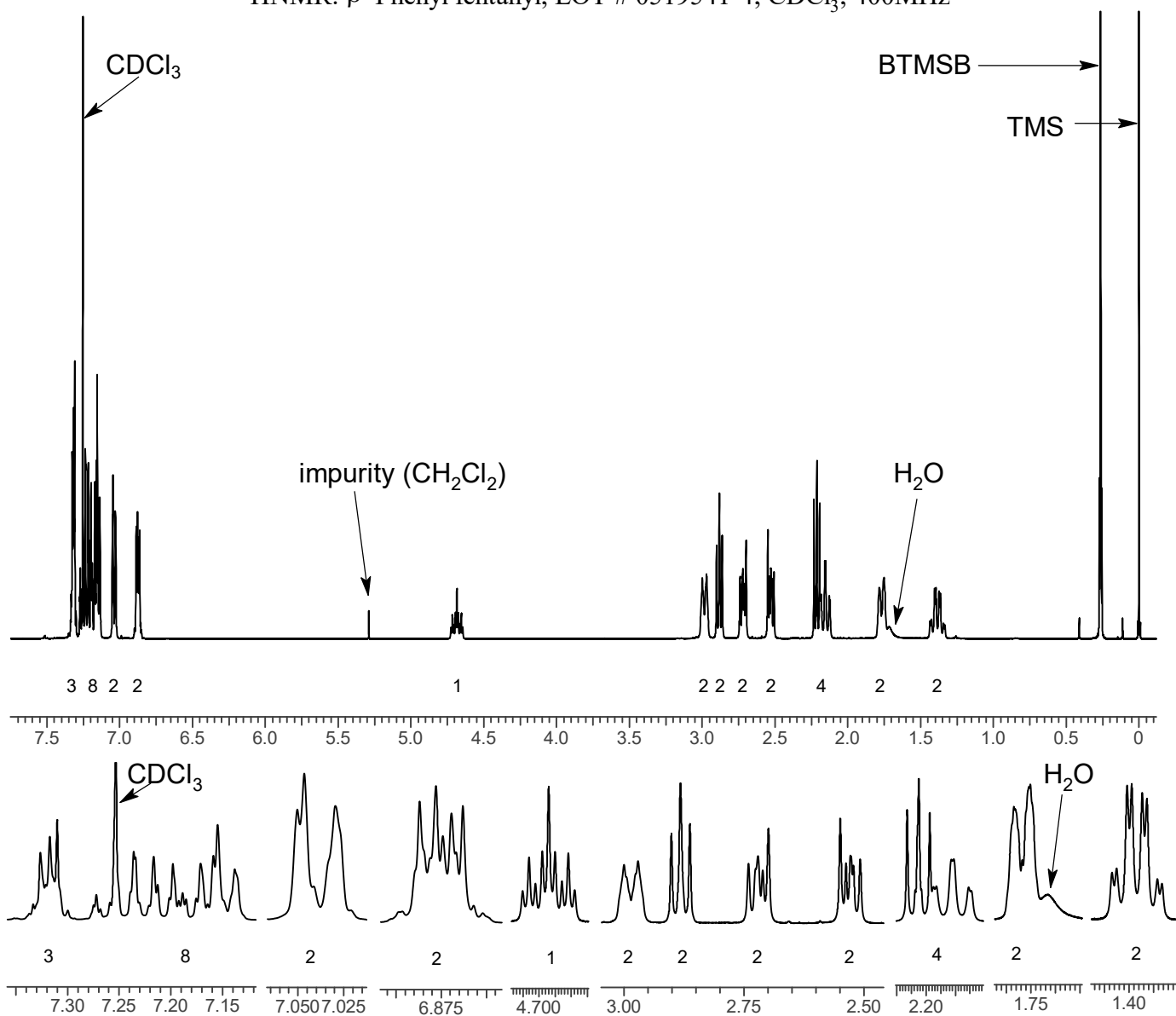



β' -Phenyl fentanyl

The Drug Enforcement Administration's Special Testing and Research Laboratory generated this monograph using structurally confirmed reference material.



3.2 GAS CHROMATOGRAPHY/MASS SPECTROMETRY

Sample Preparation: Dilute analyte ~4 mg/mL in MeOH

Instrument: Agilent gas chromatograph operated in split mode with MS detector

Column: HP-5 MS (or equivalent); 30m x 0.25 mm x 0.25 μ m

Carrier Gas: Helium at 1.5 mL/min

Temperatures: Injector: 280°C MSD transfer line: 280°C

MS Source: 230°C MS Quad: 150°C

Oven program:

1) 100°C initial temperature for 1.0 min

2) Ramp to 280°C at 12 °C/min

3) Hold final temperature for 30.0 min


Injection Parameters: Split Ratio = 25:1, 1 μ L injected

MS Parameters: Mass scan range: 30-550 amu Threshold: 150

Tune file: stune.u Acquisition mode: scan

Retention Time: 26.343 min

EI Mass Spectrum: β' -Phenyl fentanyl; LOT # 0519541-4

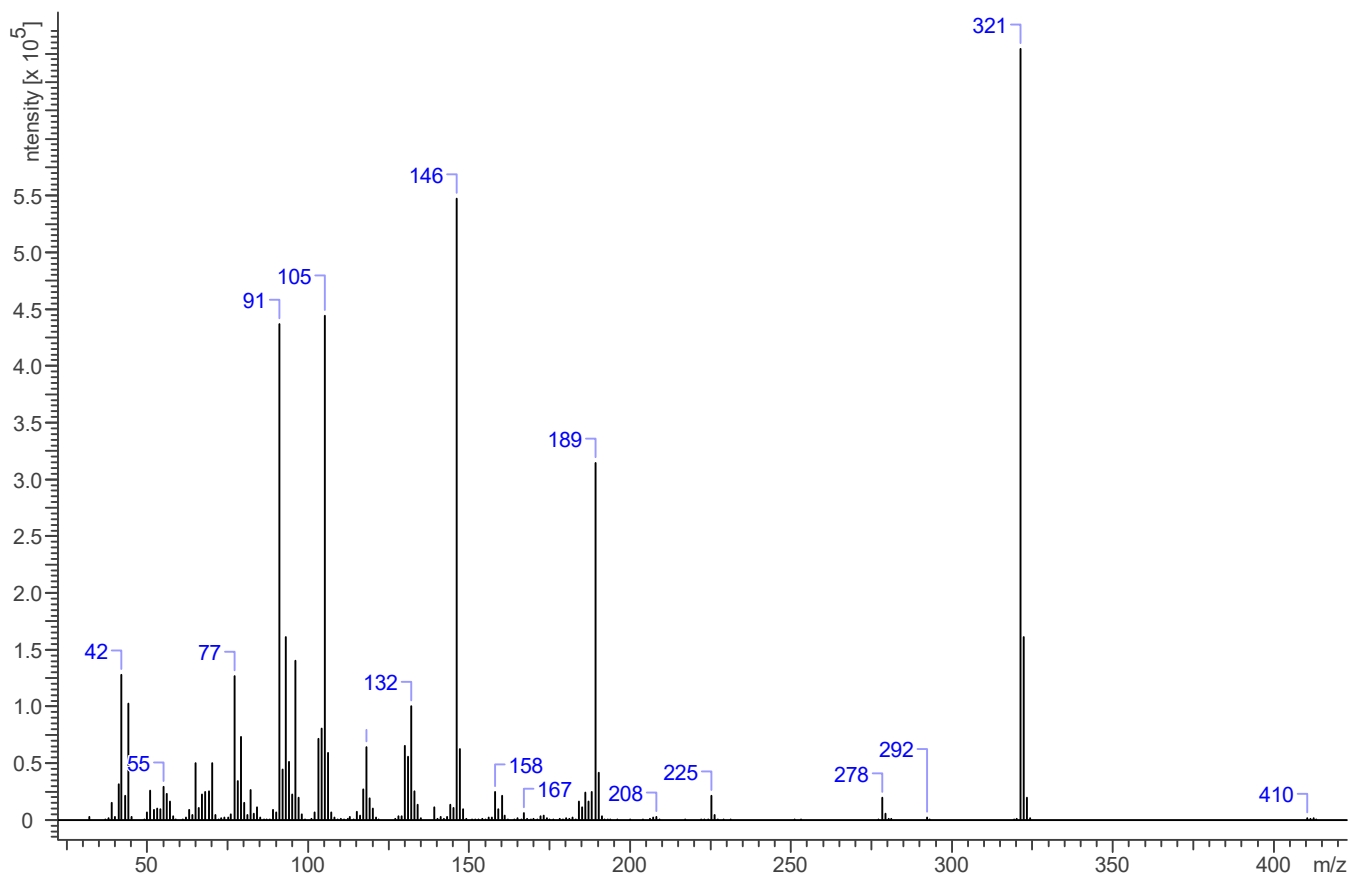



β' -Phenyl fentanyl

The Drug Enforcement Administration's Special Testing and Research Laboratory generated this monograph using structurally confirmed reference material.



3.3 INFRARED SPECTROSCOPY (FTIR)

Instrument: FTIR with diamond ATR attachment (1 bounce)

Scan Parameters:
Number of scans: 32
Number of background scans: 32
Resolution: 4 cm⁻¹
Sample gain: 1
Aperture: 150

FTIR ATR (Diamond 1 Bounce): β' -Phenyl fentanyl; LOT # 0519541-4

